


4 Types Of Waste Bins In The Laboratory

General Waste

Clinical Waste (Autoclavable Bag)

Clinical Waste (Biohazard Bag)

Sharp Bin

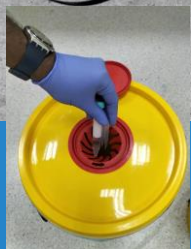

- Microorganism (Bacteria/Fungus)
- Human/Animal tissue

- Glove
- Tissue
- Disposable Pasteur Pipette

- Lancets
- Needles
- Syringes

NMF-2024

5 Steps On How To Dispose Of Biohazardous Waste In The Laboratory



1. Put the waste in an autoclavable plastic bag.

2. The waste is tied using a cable tie.

3. The waste must be autoclaved before being disposed of.



5. The waste must be put in the yellow waste bin.

4. After being autoclaved, waste must be put in a biohazard bag and tied using a cable tie.

Created By :
Norwani Mohammad
Mohd Faizul Husin
2024